



Tablet Press Controller (TPC) Short User Manual

STARTUP

You can **Login** as Operator, Supervisor, or Maintenance personnel.

The setup recipe and batch will start automatically.

To continue, adjust the press, select another recipe and start a new batch.

RECIPE

- While the batch is running, any changes by operator to recipe target or limits will go into effect **for the current side** after you click **SUBMIT** button.
- When the batch is paused, supervisor can change, save, select, or add new recipe. For a double-sided press, any changes will apply to **both press sides**.

STATUS

- User can adjust dosing in **Manual mode**.

PUNCHES

- Per punch Compression Force histogram for either side with Rejection Limits displayed.

REPORTS

- **Audit Trail Event Log** includes
 - All significant actions
 - All alarms
 - Every 15 minutes, the batch status if batch is running.
- **Batch Report** includes
 - Recipe information
 - Current batch status.

SETUP

- After “S” sequentially rejected tablets, the system will sound an alarm.
- After press starts, all tablets from “R” revolutions will be rejected.
- After “T” rejected tablets in “N” revolutions, the system will sound an alarm.
- After “M” minutes of inactivity the screen becomes frozen and requires a logon.
- After “A” revolutions in manual mode, the press will switch to control mode.

BATCH

- Batch will stop when batch size limit is reached.
- While the **batch is running**, you can
 - temporarily change recipe target and limits
 - change user
 - stop batch
 - see Audit Trail event log (Supervisor access) and report for the current batch.
- When **batch is paused**, the TPC is operating in manual mode and you can
 - resume batch
 - start a new batch
 - change batch parameters
 - change current recipe (Supervisor access)
 - save current recipe (Supervisor access)
 - select another recipe
 - create a new recipe (Supervisor access)
 - change user
 - change setup parameters (Supervisor access)
 - manage passwords (Supervisor access)

- see Audit Trail event log (Supervisor access) and report for the last batch.
- Batch **RESUME** means tablet counts are preserved.
- Batch **START** means tablet counts are reset.
- Lot number is just a label that is reported in the Batch Report.
- You can have different lot numbers for the same batch name.
- Audit Trail file is appended with new events until you start a new batch. Then a new Audit Trail file is created.

MODE OF OPERATION

The mode button in the header will toggle between Manual and Control mode.

- In Manual Mode:
 - The dosing control is disengaged
 - The tablet rejection is disabled
 - Automatic shutdown is disabled
 - User can start/stop press
 - User can adjust dosing cam from the **STATUS** screen

After “A” revolutions from the batch start, the system is automatically switched to control mode. The number “A” can be changed from the **SETUP** screen.
- In Control Mode:
 - The dosing control is engaged
 - The tablet rejection is enabled
 - Automatic shutdown is enabled. Press will shut down when batch limit is reached, on a critical alarm, when average force is outside shut-down limits, when individual punch force is outside the limit
 - User cannot adjust dosing cam from the **STATUS** screen.

PASSWORDS

- Operator can
 - Select recipe
 - Change batch name, size, lot number
 - Start / stop / resume batch
 - Toggle manual and control modes
 - Change recipe target and limits on the fly
 - Generate Batch Report
 - Adjust dosing cam (in manual mode)
 - Adjust press speed
 - Copy log and report files to diskette
 - Shut down the computer
 - Logout by clicking on the **USER** bar on the header.
- Supervisor can
 - Do everything an operator can do
 - Change, save, select, and create new recipe
 - Change, save, and create new users and passwords
 - Modify setup parameters
 - Exit the program
- Maintenance user can
 - Do everything a supervisor can do
 - Modify calibration factors